



Valiteq Glove Fingertip & Media Fill Test Kit Instructions

The AAOA offers resources and tools to help you meet the USP General Chapter <797> compliance criteria for sterile compounding.

The [USP General Chapter <797> Sterile Allergen Extract Compounding Online Module](#) offers you and your practice a resource to meet the compounding personnel criteria. USP General Chapter <797> requires personnel to demonstrate proficiency in knowledge of principles and skills for sterile compounding every 12 months and to document training and competency. Using an adaptive learning platform, the AAOA's online module is a great tool to help assure you can meet and document the proficiency defined in USP <797>. As an adaptive learning tool, participants will succeed through the learning process. It is a great to both learn about sterile compounding, as well as document proficiency.

AAOA has also secured an AAOA members discount with [Valiteq](#) for its testing kits (AllerTEQ AL1) used to meet the media fill and gloved fingertip testing requirements. AAOA members will get a 15% discount on the kits. Each test kit requires a 2-week incubation process, 1-week growth promotion, and a written report at 4 weeks. To order your test kits, please fill out an online [Order Form](#) linked below and use the code AAOAUSP2020. Email the Order Form to aaoa@valiteq.com. Invoices are issued upon test kit delivery. Valiteq has additional test kit options and tools that go beyond the USP requirements to help ensure staff understand sterile preparation. Feel free to enquire.

AllerTEQ AL1

- Initial test set (AL1DI) **\$187 AAOA member pricing** plus shipping (\$220 list price plus shipping for non-AAOA members)
- Initial test set (AL1DI) includes three sets of (2) contact plates for the triplicate glove fingertip garbing test and one set of (2) contact plates for the post media fill glove fingertip testing as well as the media fill test
- Requalification test set (AL1DQ) **\$172.55 AAOA member pricing** plus shipping (\$203 list price plus shipping for non-AAOA members)
- Requalification test set (AL1DQ) includes two sets of (2) contact plates, one set for the glove fingertip testing and one set for the post media fill glove fingertip testing as well as the media fill test kit.
- [AllerTEQ AL1 Order Form link](#)


Valiteq Glove Fingertip & Media Fill Test Kit Instructions

Receiving and handling your test kits

Your shipment arrives by second day air and it contains perishable contents, so it's important to follow the instructions on how to unpack, store and return the test kit.

Notice - Do not discard any of the contents, until the test is completed and ready to be returned Valiteq for analysis.

1. Content sheet

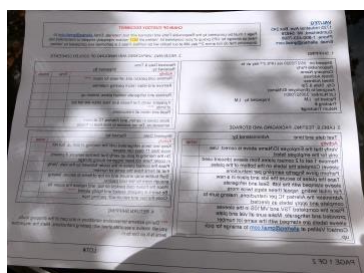

- Describes what is in the test
- Denotes employee doing the test
 - Initial each activity

This form needs to be returned along with the test once the test is completed.

2. Unpacking Instructions




Includes unpacking instructions for normal weather conditions as well as for warm weather shipping & packaging (plates need to be refrigerated). Please read the 15 steps before unpacking and storing your test kit to ensure unpacking and storing is done correctly. Also, you'll find the return instructions.

3. Test kit and plates:



Your shipment contains two boxes. The smaller one is the Media fill test kit. The larger box labeled Perishable contains the 3 plates (follow instructions on how to correctly store).

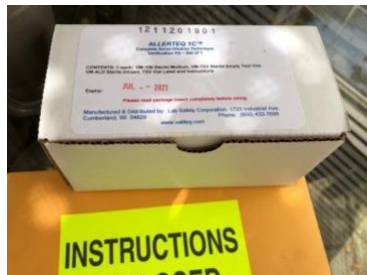

Media fill test kit (vials, instructions and shipping information)



3 Testing Plates – Need to be refrigerated

If you have any questions, please check out our [Unboxing Video](#) or call our office at 202.955.5010.